

Demo - How to use the MicroCT analysis computer

By PCMD MicroCT Imaging Core


Youtube link: <https://www.youtube.com/watch?v=qHHcB6KJJe4>

All our video tutorials are listed on our official website


<https://www.med.upenn.edu/orl/uct/>

Before you begin, we strongly suggest you restart this computer!


Click the **Start** button > select "Restart" in the menu.

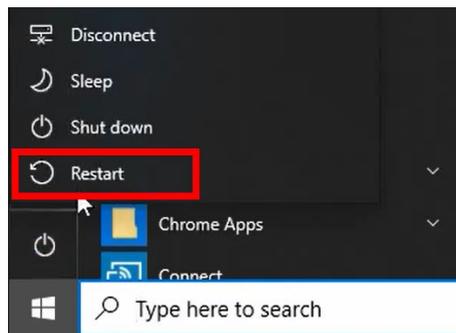

Important! Please make sure **Caps Lock**, **Num Lock**, and **Scroll Lock** are all OFF on the keyboard!

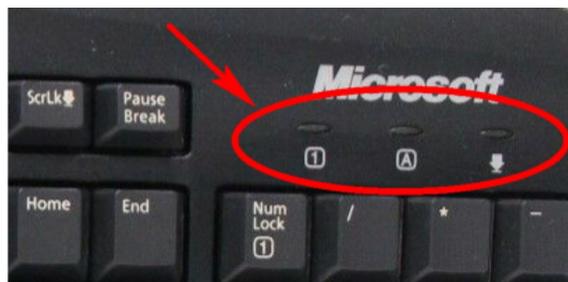

Suppose you want to do analysis on MicroCT35. Double-click the icon "uCT35":

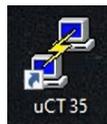

A command window will pop up (black background).

Note: If you want to analyze samples in other scanners, make sure you are accessing the correct scanner for analysis: uCT45, vivaCT40, vivaCT80

```

demo.yiluzhou.com - PuTTY
Using username "microct".
Pre-authentication banner message from server:
| Welcome to HP OpenVMS Industry Standard 64 Operating System, Version V8.4
|
| End of banner message from server
|
| Last interactive login on Monday, 23-MAR-2020 18:40:06.31
| Last non-interactive login on Wednesday, 25-MAR-2020 00:38:16.17
DISK1:[MICROCT]
$

```

For evaluation analysis, type `uct_evaluation` in the command window then press `Enter`. The evaluation window will pop up in a few seconds.

```

demo.yiluzhou.com - PuTTY
Using username "microct".
Pre-authentication banner message from server:
| Welcome to HP OpenVMS Industry Standard 64 Operating System, Version V8.4
|
| End of banner message from server
|
| Last interactive login on Monday, 23-MAR-2020 18:40:06.31
| Last non-interactive login on Wednesday, 25-MAR-2020 00:38:16.17
DISK1:[MICROCT]
$ uct_evaluation

```

If you want to open another evaluation window at the same time (e.g.: you need to compare images), you may double click "`uCT35`" to open a second command window.


For microCT 3D images, type `uct_3d` in the command window, press `Enter`.

```

demo.yiluzhou.com - PuTTY
Using username "microct".
Pre-authentication banner message from server:
| Welcome to HP OpenVMS Industry Standard 64 Operating System, Version V8.4
|
| End of banner message from server
|
| Last interactive login on Monday, 23-MAR-2020 18:40:06.31
| Last non-interactive login on Wednesday, 25-MAR-2020 00:38:16.17
DISK1:[MICROCT]
$ uct_3d

```


If you encounter software problems, please **simply restart this computer** or close all command windows (black background). Then, click taskbar options, right click on the icon labeled "Xming: Server:0.0".

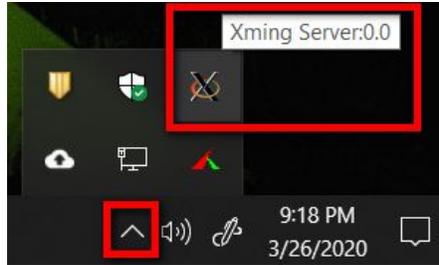

Select "exit".



Select "exit" again in the prompt window.

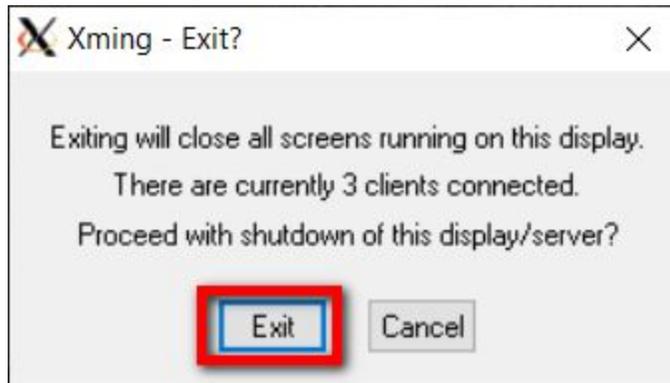

Double click the "MicroCT Preload" on the desktop:



Then you can resume "uCT35" for analysis (step 1).

If issues still persists, please take screenshot and include a brief description of the issue(s) to pcmd.microct@gmail.com